

LIMITED WARRANTY

The manufacturer guarantees that this unit is free from defect in materials and workmanship when it leaves the factory and undertakes to replace or repair the unit if it proves defective in normal use or during servicing for a period of 1(one) year. The date of original installation and is for the benefit of the original purchaser only. The liability under this warranty is limited to repairing the defective unit or any part of the unit provided it is sent carriage paid to an authorized dealer. All other Warranties ,expressed or implied , statutory or otherwise, including without limitations any implied Warranty of Merchantability or fitness for purpose are excluded .The Seller shall in no event be liable for direct,indirect or consequential damages in connection with the products.

This unit is at all times to be used according to the instruction manual and for its normal purpose.

This Warranty is not effective if damage occurs because of accident, carelessness,improper installation,lack of proper set-up,supervision when required or if the equipment is installed or operated in any manner contrary to the installation and operating instructions. In these cases,repairs will be made at a reasonable cost. Work performed by unauthorized personnel or unauthorized service agencies voids this Warranty.



INSTRUCTION MANUAL MODEL 723 MIXER



ELECTRICAL CONNECTIONS

1. The appliance operates on 220-230V/50Hz direct current. It must be plugged into a safety socket
2. The technical characteristics of the appliance are given below:

TYPE:	723
POWER:	350 Watt
MAINS VOLTAGE:	220/230V - 50Hz
WEIGHT:	6.14 kg
HEIGHT:	0.46 m
CATEGORY:	TYPE 1
FIRE PROTECTION RATING:	TYPE B

ATTENTION

1. Read through the instructions carefully before trying to operate the appliance.
2. Keep the instructions in a safe place for future reference.

SAFETY INSTRUCTIONS

1. Do not cut off the plug, because the appliance is not designed for permanent electrical connection.
2. Do not use an adapter or an extension cord to connect the appliance.
3. The lead must not be allowed to touch hot surfaces or sharp corners
4. Do not try to run the appliance if it should fail to operate properly or if the lead appears to be defective, for this may cause fire or electrical shock. ONLY authorised service agents are permitted to repair or change parts of the appliance.
5. never move the appliance by pulling on the lead
6. Do not run the appliance if your hands are wet.
7. **Do not submerge the appliance in water or any other liquid.**
8. Do not touch the rotating parts of the appliance as this may cause minor injuries.
9. Be sure to keep the appliance clean. The rotor and container are made of stainless steel. Since these are the parts of the appliance that come in direct contact with food, they must be kept scrupulously clean.
10. Never try to clean the appliance while it is plugged in.
11. Do not allow small children or handicapped persons to operate the appliance.
12. Do not disassemble the appliance.

BEFORE USING FOR THE FIRST TIME

1. Unwrap the appliance and wash the container, using an ordinary washing-up liquid and being very careful to rinse thoroughly.
2. Wash the rotor by rubbing with a soft soapy cloth. Rinse with a damp cloth.

OPERATING INSTRUCTIONS

1. Place the appliance on a dry, stable, level surface.
2. Plug the appliance in.
3. Place ingredients in the special container. The appliance is designed for use with liquids only. Do not attempt to use this appliance to crush ice or shred other solids.
4. Set the switch on the head of the appliance to the desired rotor speed (0-I-II). Insert the rotor shaft into the container, set the container on its base and press down to establish contact. There are three positions available to accommodate 11 - 17.5-cm glasses. Select one of the positions by pressing down. The appliance will continue to run as long as you continue to press down.
5. It is recommended that you wash the appliance after each use.
6. You should unplug the appliance if it is not going to be used for a considerable period of time.

CLEANING THE MIXER

1. Unplug the appliance
2. Wash the container, using an ordinary washing-up liquid. This will ensure compliance with all health regulations.
3. Clean rotor and other stainless steel parts by rubbing with a soft soapy cloth. Rinse with a damp cloth.